

January 2019

*Preparing Samples of Cannabis Plant for THC/CBD Analysis*

Porvair Sciences reports on a highly effective protocol, using its Plant Genomics microplate, for preparing cannabis samples for analysis of Tetrahydrocannabinol (THC) and Cannabidiol (CBD).

With increasing legalization of cannabis for medicinal and research purposes, demand for accurate and high throughput analysis of its active ingredients has dramatically risen. Cannabis plant material contains several cannabinoids for which the human brain has evolved receptors. The two most important are the psychoactive THC and the analgesic CBD which has therapeutic uses. The ratio of THC and CBD is crucial for differentiating medicinal versus recreational cannabis. This ratio is important for taxation authorities, medical regulators and growers alike.

Accurate determination of the THC/CBD ratio is critically determined by effective sample preparation of the samples of cannabis leaf, bud and flower. Traditional sample preparation methods have used bead-beating or grinding mills to homogenise these materials. When conducted in polypropylene microplates the extreme forces applied by these machines often cause damage to the plates leading to cracking, leaking and ultimately, cross-contamination.

With more than thirty years' experience in microplate manufacture, Porvair Sciences has applied modern design concepts and mould analysis to this problem to create a solid, stable and strong microplate that can withstand the applied forces in bead-beaters whilst maintaining a standard ANSI/SLAS footprint and key dimensions. By adding reinforced walls to the bottom of the well and backing these with supporting ribs on the underside of the plate, Porvair Sciences has produced a 2 ml square well block (Plant Genomics plate) that can be used safely and with confidence for Cannabis plant extraction. The Plant Genomics plate has a generous working volume of 1.90 ml per well in a 45mm high plate. The polypropylene used to mould the plate is pure, virgin, extractable free polymer, rigorously tested for leachates thereby ensuring no contamination from the plate will affect your THC/CBD analysis.

In the described protocol, sample preparation using the Plant Genomics plate in a GenoGrinder takes just 2 minutes. Following this, the dry ground Cannabis powder is transferred into the wells of a Porvair P3 filter microplate for extraction under agitation with methanol or isopropanol to dissolve the oily residues in the plant material. After agitation, the P3 plate is placed onto a Porvair Microlute manifold and vacuum applied to draw the purified extract down through the filter and into the wells of a collection plate. The collection plate containing the organic extract is then taken for HPLC analysis to determine the THC/CBD ratio. To determine total THC and CBD in the Cannabis plant sample, the purified extract should be gently evaporated to dryness using a Porvair MiniVap blowdown evaporator. The sample should then be redissolved in a known volume of solvent ready for HPLC analysis. Both methodologies are reported to be straightforward and repeatable.

For a copy of this report and further information on the Plant Genomics plate please visit <https://www.microplates.com/resources/porvair-seed-genomics-plates-flyer/> or contact Porvair Sciences Ltd on +44-1978-666222 / +1-800-552-3696 / [int.sales@porvair-sciences.com](mailto:int.sales@porvair-sciences.com).

Established in 1992, Porvair Sciences is one of the largest global manufacturers of ultra-clean microplates and top-quality microplate equipment for life science, synthetic chemistry and many other applications. Porvair Sciences Ltd. is a wholly owned subsidiary of Porvair plc.

-----



**Worldwide HQ**

**Porvair Sciences Ltd**

Clywedog Rd South  
Wrexham Industrial Estate  
Wales LL13 9XS  
UK

Tel: +44 1978 666222

Email: [int.sales@porvair-sciences.com](mailto:int.sales@porvair-sciences.com)

Web [www.porvair-sciences.com](http://www.porvair-sciences.com)